

Working safely at safety cabinets

1



Observe user information!

2



Instruct and authorize personnel!

3



Safety cabinet:
Single workplace!

4



Use personal protective
equipment!

5



Safe: Green control lamp!

6



Unsafe: Red control lamp
+ audible warning signal!

7



Critical activities only with front
screen in working position!

8



Do not block air intake
openings!

9



Do not block exhaust air
opening!

10



Safe area: closed work
surface!

11



Workspace: No storage
area!

12



No strong heat sources (e.g. gas
burner) in the workspace!

13



Organise transport of 12
material well!

14



Clean and disinfect material
before taking in or out!

15



Close communication between
pharmacy and station!

16



Quality-oriented production!

17



Clean and disinfect workspace
daily!

18



Keep to good microbiological
and/or aseptic techniques!

19



Avoid quick arm movements!

20



Avoid disturbing the air flow in
the surrounding area!

21



Generally only particles
but no gases are deposited!

22



Regular inspection by
authorised and competent
service personnel!



BERNER INTERNATIONAL GMBH
Mühlenkamp 6 • D-25337 Elmshorn
Germany
www.berner-international.eu
info@berner-international.de
T.: +49 (0) 41 21 / 43 56 - 0
F.: +49 (0) 41 21 / 43 56 - 20

BERNER
safety systems
made in Germany

More information: www.berner-international.de/safety_cabinets